

Workflow Addon API

The WF API works on the same logic as the LabCollector API. Please see our [knowledgebase on LabCollector API](#).

Required headers:

- X-LC-APP-Auth (authentication)
- Accept (return type)

PUT Actions:

execute_job

Used to start an job in the WF addon. This start has all the same functionalities as a normal job start from the WF addon (like automatic start).

Required Parameters:

1. name (name of new job)
2. template (template of new job) -> can be the template name or the template ID

Example:

```
http://[my_labcollector_url]/extra_modules/workflow/webservice/?action=execute_job&name=XXXX&template=XXXX
```

complete_node

Used to trigger the completion of node API Execution type. Or instead setting the node to error status.

Complete:

Required Parameters:

1. job (name of job where the node is) -> can be the job name or the job ID
2. node_id (node identifier) -> always a negative number (can be seen in the API Execution node in the template side)

Example:

```
http://[my_labcollector_url]/extra_modules/workflow/webservice/?action=complete_node&job=XXXX&node_id=-XX
```

Error Status:

Required Parameters:

1. job (name of job where the node is) -> can be the job name or the job ID
2. node_id (node identifier) -> always a negative number (can be seen in the API Execution node in the template side)
3. obs (when setting a error status observations related to that error are mandatory)

Example:

```
http://[my_labcollector_url]/extra_modules/workflow/webservice/?action=complete_node&job=XXXX&node_id=-XX&status=E&obs=XXX
```